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August 2025

Photometry designed as a system – why it is not recommended to use third-party reagents on Hach instruments

System calibration to eliminate errors

Hach reagents and instruments are developed and designed to work as an integrated system, eliminating calibration and variability errors from the outset to obtain best possible results.

Maintaining quality by avoiding negative impact of raw materials

It is not just a one-time effort to calibrate instruments with reagents. Hach reagents are produced daily, and it is an ongoing process to always maintain the quality and analytical performance of the photometric system. In an ever-changing chemistry raw material market, variation in raw materials may have a negative impact on quality of results. Thus, stringent internal quality control checks are required to ensure that each single production lot of each reagent matches with the original Hach calibration curves within given tolerances. Hach's quality management system ensures that Hach reagents and Hach instruments work best together – every day. And in case raw material variations are significant and a new calibration curve is required, the Hach R&D team takes care of this quickly. The new calibration data and lot-numbers of updated reagents will be communicated to Hach users to ensure that instruments and updated reagents continue to match correctly. This Hach service to maintain best possible results is not feasible with third-party reagents.

Hach uses special reagent compositions to increase stability and to reduce interferences

Besides common or mandatory components for the photometric color reaction, Hach formulations often contain special additives. These enable long product life and often help receiving best possible results after shipping and storage. In addition, such additives may reduce the effects of common sample interferences. It is unknown if third-party reagent suppliers always use the same reagents composition as Hach. Deviations in formulations may have an impact on reagents shelf life or may lead to different results in specific sample matrices.

Hach analytical support is not feasible for third-party reagents

In case analytical support is required, the Hach technical support team usually asks for the following information to provide help: "What instrument, which Hach reagent and which lot number did you use". This holistic information about the photometric system helps to identify the root cause and to solve the problem. Such analytical support can hardly be provided in the case of third-party reagents being used instead of the entire Hach system. It is then difficult to identify the root cause of a problem and whether the instrument or the reagents do not work correctly. Hach disclaims all liability and provides no warranty whatsoever with respect to such third-party reagents.

CARSTEN SCHULZ

Global Product Manager

Hach Company | www.hach.com